

## Instructions for Uploading Firmware and Software

Attached are the 2 files for the miniGC for software and firmware.

To install the software. Close the software and run the .exe file.

To install the firmware. FIRST download the .bin file the Local Disk C drive. There you will see a previous version of the .bin file. Add the new .bin file there.

Turn off the miniGC.

Attach the USBmicro cable between the back of the system and the laptop. The micro end of the cable goes into the labelled port on the back of the system (miniGC or miniLC) and the USB side of the cable connects to a USB port on the laptop.

Navigate in software

- 1 click on the menu Icon
- 2 Click on Instrument
- 3 Scoll down to GC Firmware
- 4 Click update and find the file in the C drive.

Once finished unplug the USB cable from the laptop or GC before powering on the miniGC.